

Paraffin dispenser

User manual



Attention: Before using the machine, please read this Instruction and keep the Instruction for further reference.

Warning

To prevent fire or shock hazard, do not expose this unit to rain or moisture. This equipment should be used with 110V, 60Hz.

No user-serviceable section inside. Refer servicing to qualified service personnel.

Application


This paraffin dispenser is designed for melting paraffin and dispensing the paraffin. The sole purpose of this paraffin dispenser is for histology use, research, teaching purpose. The Paraffin Dispenser is not recommended for any other uses than the ones described below in this user manual.

Assembly Instructions

1. Remove Paraffin Dispenser from styrofoam packing and place it on a stable worktable. Remove all plastic bags and paper covering (these can be discarded).
2. Connect the Paraffin Dispenser power supply to an 110V outlet and your Paraffin Dispenser is ready for use.

Operation

1. Prior to use, fill the Paraffin Dispenser with solid paraffin at least half-full. It's also recommended not fill the Paraffin Dispenser with too much paraffin for it might overflow. HAZARD-Having liquids near or in contact with power supply may cause electric shock.
2. After making sure that the Paraffin Dispenser is plugged in, turn on the machine with the switch in the back. The switch indicator light (#3) should light up to indicate the machine is working.
3. Before using, check the temperature to make sure that your set temperature is above the melting point of your paraffin. (Your paraffin box should indicate melting temper).
4. To adjust the temperature, press the set button (#4), then use the up and down arrow button (#5) to set the temperature at the desired setting. This adjustment should be visible in the setting display panel (#2). The actual temperature of the wax basin can be seen in the actual temperature display panel (#1). Please wait patiently for the Paraffin Dispenser to melt the wax before it starts dispensing.
5. Once the Paraffin Dispenser reaches your set temperature, the heat light indicator (#3) will go out.
6. Timer function. You can set the heating time with the timer. Push the set button (#4) and the time can be set on the setting display (#2). Then use the up and down arrow buttons (#5) to set the timer. This adjustment is visible in the setting display panel (#2). Once you reach the desired time, hit the set button (#4). Your timer will count down and buzz (meaning timer is off).
7. The faucet heating intensity can be adjusted with the knob at the rear of the device.

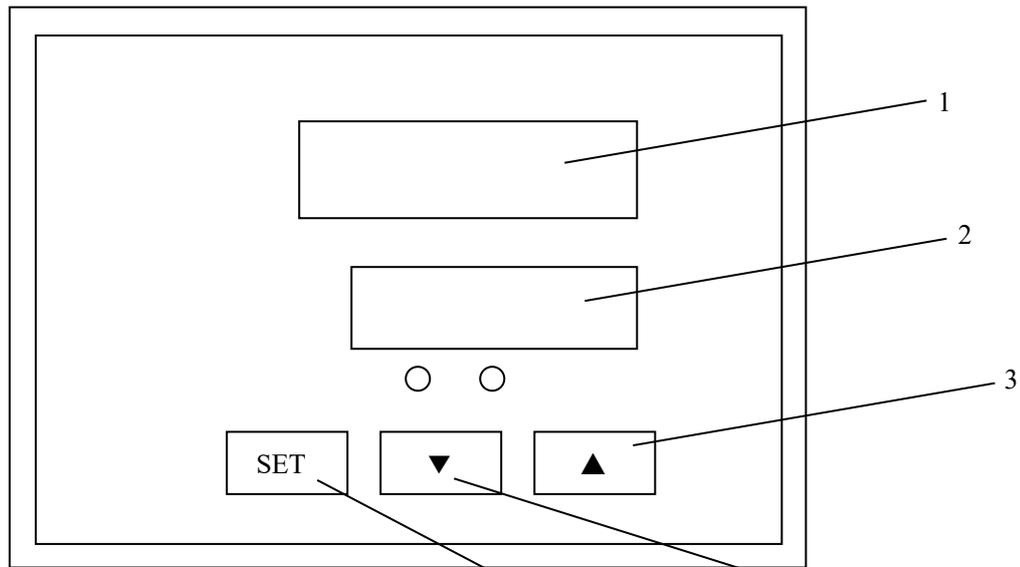

- 1) Temperature Display
- 2) Setting Display
- 3) Switch Light indicator
- 4) Set button
- 5) Up and Down arrow Buttons

Cautions

1. Unplug the Paraffin Dispenser when not in use.
2. Do not use the faucet when the temperature is below the melting point of the paraffin.
3. Be very careful with the faucet, as it may be hot because it can absorb the heat from the Paraffin Dispenser when the dispenser is set high temperature.
4. Please contact your dealer with any operating problems. Only authorized repair technicians should perform repairs.

Maintenance

1. Always unplug Paraffin Dispenser before moving or cleaning. HAZARD-Having water or any other liquid near or in contact with power supply may cause electrical shock.
2. Empty and clean Paraffin Dispenser regularly. To clean use a lint free type of towel and sponge when necessary. Avoid using any type of instrument/tool/chemical that can scratch or damaged the Paraffin Dispenser.
3. The Paraffin Dispenser has been tested and inspected before shipment.